

Muffle Furnace

User manual

SAFETY INFORMATION

1. Heating ramp rate should not exceed 5°/min.
2. Remember to wear heat protecting gloves.

Don't use without proper training.

Operation:-

1. Ensure that the area is clean and the calibration due data of the temp. Indicator is within calibration period.
2. Ensure that the instrument is connected to the power supply.
3. Switch on the main power supply, glowing of red lights at mains indicates the power supply.
4. Switch on the instrument by ON position which leads to activation of green control bulb and the temp controller.
5. Set the temp required by pressing and holding v buttons and then ENT key is pressed to finish operation.
6. The digital display shows the actual temp of furnace and the red light of temp controller glows
7. When the temp reaches at the setting point, the red light of temp controller switched OFF and green light will glow
8. The equipment is now ready for operation.

