

Programmable oven

User manual

SAFETY INFORMATION - The programmable oven is designed for safe and efficient operation when used properly and in accordance with this manual. Failure to observe the following precautions could result in serious personal injury:

- ❖ The programmable oven is an electrical instrument. To avoid electric shock, please observe all standard precautions.
- ❖ All the connections between different parts of instrument should be proper.
- ❖ **Don't use without proper training.**

Operation:-

1. Check all connections carefully.
2. Switch on main power.
3. After few seconds, the current temperature inside the chamber. will display on the PID controller screen.



4. We can set our programme in maximum 8 steps in PID controller. Kindly make the programme accordingly.
5. For enter the programme, press the SET button 7 times and screen will display as following image:



6. From here, enter the programme steps in PID controller.
7. After entering the program steps, come back on the home screen and press the upper arrow key as shown in the image.



8. Now the programme will start. Switch on the fans.
9. Don't disturb the programme and chamber while programme is running.